

Date of Analysis: 30 September 2009

Peak Identity	Retention Time (min)	Purity (% Total Area) <sup>a</sup>
Oxadiazon	23.24	100.00

**GC/MS Instrument Parameters:**

Instrument / Ionization	Gas Chromatograph with Mass Spectrometer / Electron Impact
Solvent	Acetone
Column	J&W Scientific HP-5MS, 30 m x 0.25 mm ID, 0.25-µm film thickness
Carrier Gas	Helium at 1.0 mL/min
Oven Program	35°C, hold 2 min; ramp @ 10°C/min to 310°C, hold 7 min
Source Temperature	230°C
Auxiliary Temperature	150°C
Scan Range	25 – 500 amu
Injector Temperature	250°C
Injection Volume / Mode	1 µL / Split (100:1)
Data Analysis Software	MSD Chemstation, v D.03.00.SP1 and NIST Library v 2.0f, build 10/8/2008

